


PREPARATION

Adventurer Pro balances have two calibration methods: Span and Linearity. Some models have Internal Calibration (InCal™) for Span calibration.

- ✓ Ensure that the scale is level, stable, and undisturbed during calibration.
(Otherwise the process could fail.)
- ✓ Allow scale to warm up for five minutes after stabilizing to room temperature.
- ✓ Use suitable calibration masses (available from Ohaus).
Exact weight value depends on model. Follow prompts for indicated values to use.
(*Weight values here are examples only.*)
- ✓ Calibration may be locked out. To unlock it, use the Unlock menu. If the scale is set to Legal for Trade operation, a seal must be broken to unlock it. (*See the Instruction Manual.*)


Note: To cancel calibration, press:

INTERNAL CALIBRATION *(No external weight needed.)*

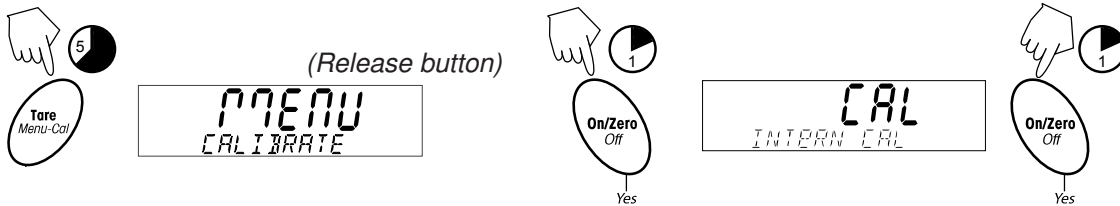


When calibration finishes, balance returns to currently selected mode.

SPAN CALIBRATION

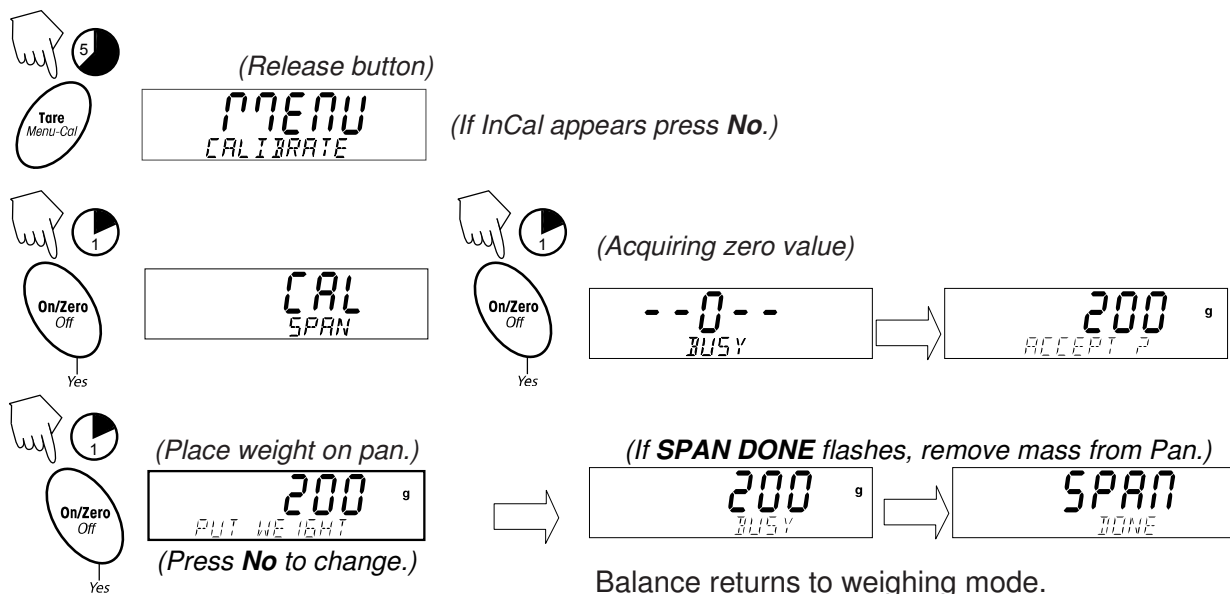

LINEARITY CALIBRATION:

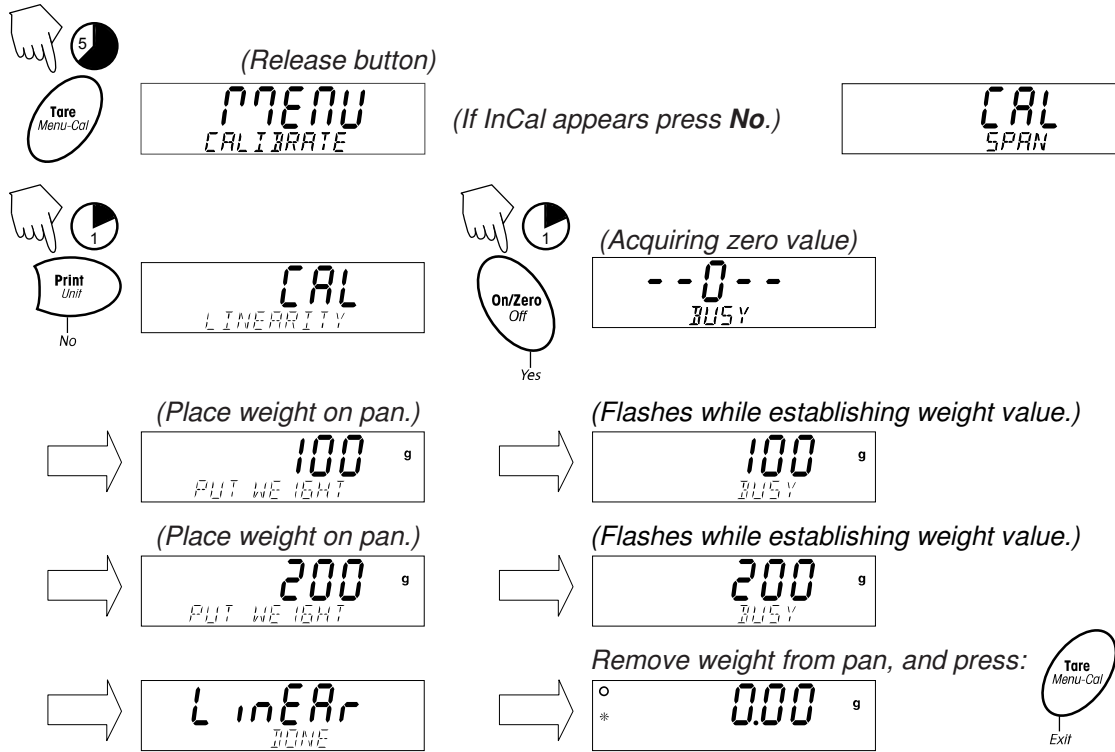

P/N 80253055