


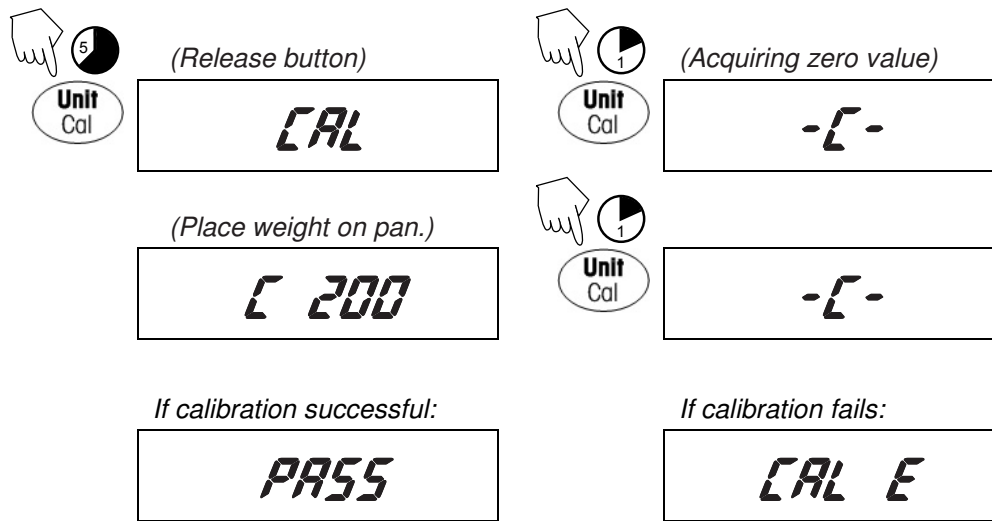
## PREPARATION

CS Scales can be calibrated as follows. Follow the prompts for calibration mass values.

- ✓ Ensure that the scale is level, stable, and undisturbed during calibration.  
(Otherwise the process could fail.)
- ✓ Allow scale to warm up for five minutes after stabilizing to room temperature.
- ✓ Use suitable calibration masses (available from Ohaus).

**Note:** To cancel calibration, long-press: 

## CALIBRATION



If **CAL E** appears, turn the scale off and try again.  
Be sure to have the appropriate calibration weight.



P/N 80253053