


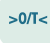


PREPARATION

Pioneer balances have two calibration methods: Span and Linearity. Some models have Internal Calibration (InCal™) for Span calibration.


- ✓ Ensure that the scale is level, stable, and undisturbed during calibration.
(Otherwise the process could fail.)
- ✓ Allow scale to warm up for five minutes after stabilizing to room temperature.
- ✓ Use suitable calibration masses (available from Ohaus).
Exact weight value depends on model. Follow prompts for indicated values to use.
(*Weight values here are examples only.*)
- ✓ Calibration may be locked. To unlock it, use the Unlock menu. (*See Instruction Manual.*)

Note: To cancel calibration, press:    until scale turns off.


INTERNAL CALIBRATION *(No external weight needed.)*

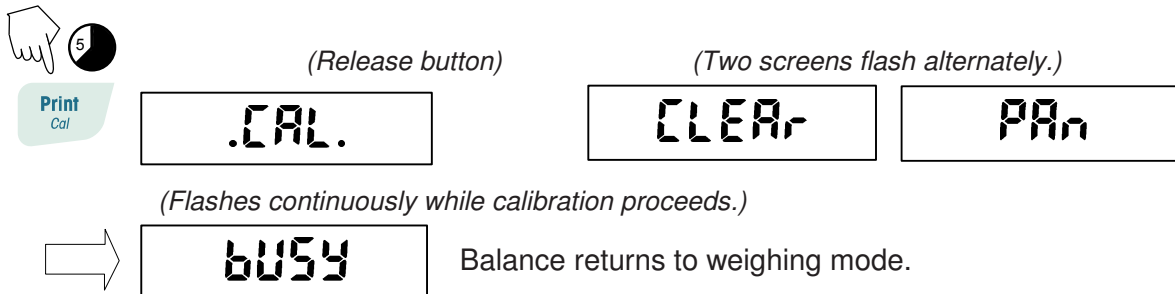

SPAN CALIBRATION

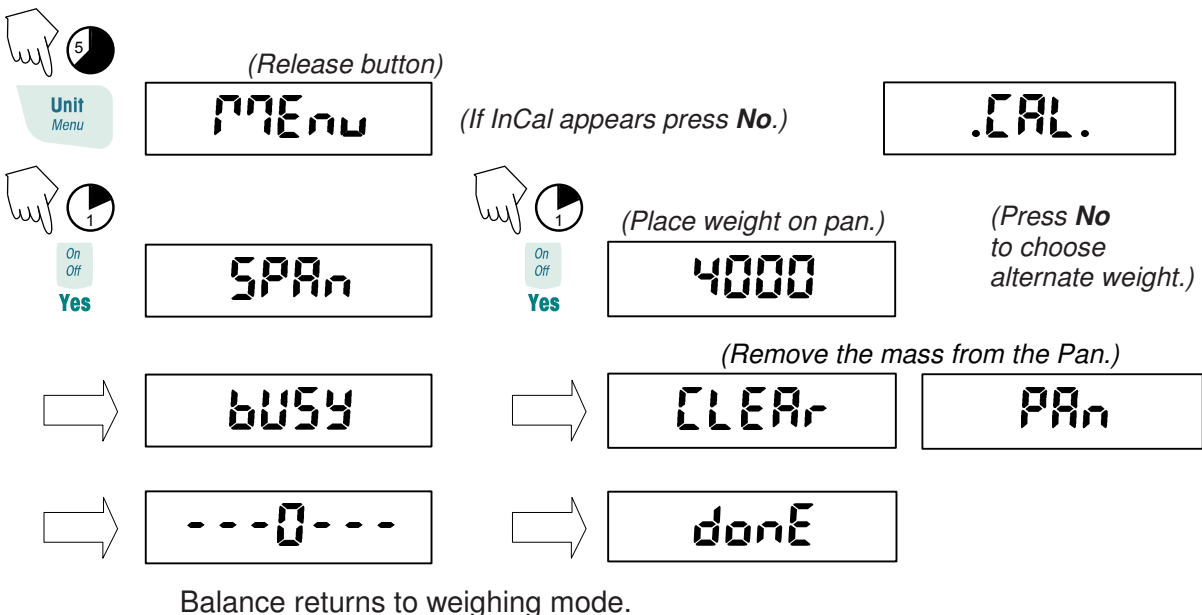

LINEARITY CALIBRATION: *(not available on InCal models)*

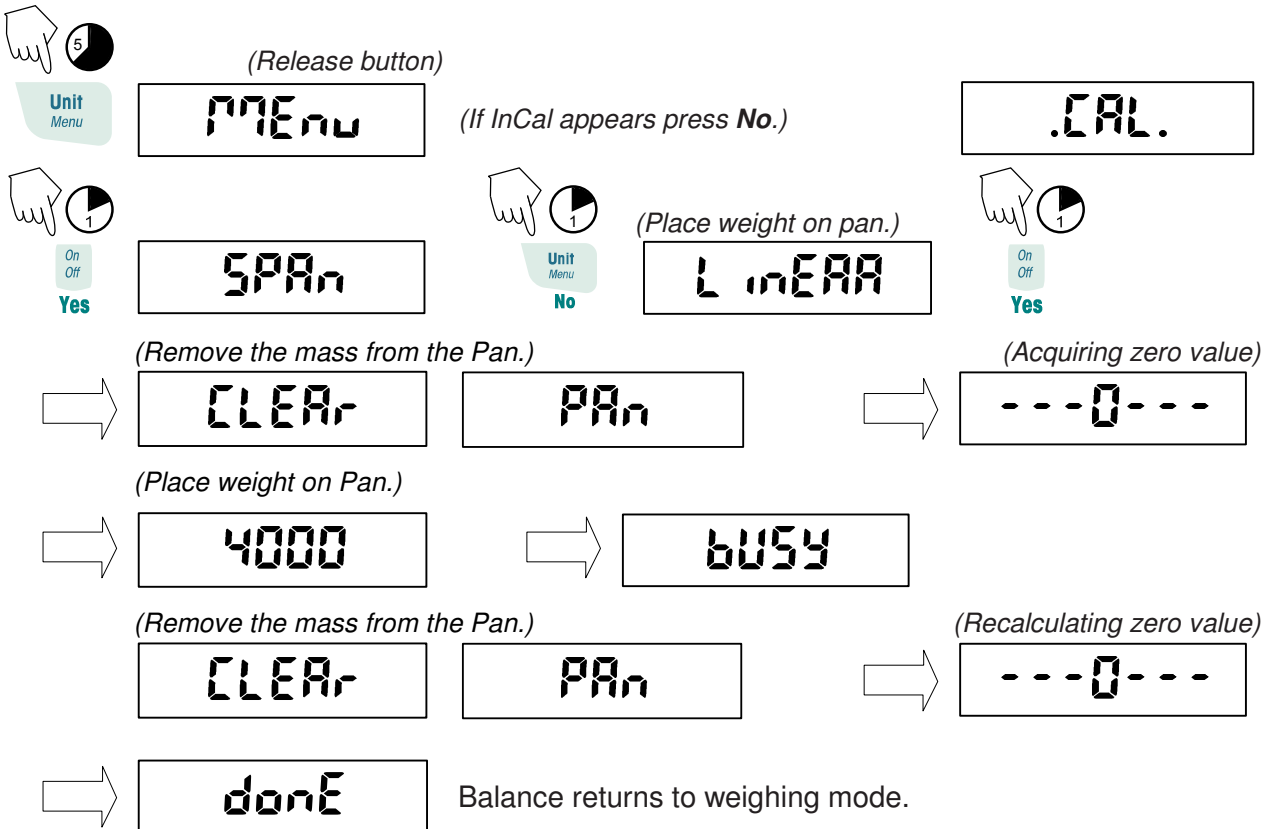

P/N 80252974